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Manual of High Speed Centrifuge Machine (Model: SRX-201)

Attention:

- Difference of weight of two diagonally positioned tubes should not be more than **0.2g (very important)**.
- Don't damage the rotary part and the central bar **(very important)**.
- Wipe out the inside water droplets after use.
- Carry out the experiment with your desired temperature but can NOT change the machine's default temperature which is **4.0 °C**.
- After finishing everything turn off the machine and also turn off the main switch from the board.

Please write note in the note book before and after use.

Keep the machine good by proper handling or don't break it by using it in the wrong way.

How to use the machine:


- (1) Turn on the main board switch then machine switch and open the cover.

- (2) Take the suspension in two centrifuge tubes then cover the tubes and balance as shown below:

- (3) Check the rotary part by small rotating by hand and check the number for this rotation on the display of machine ie.211.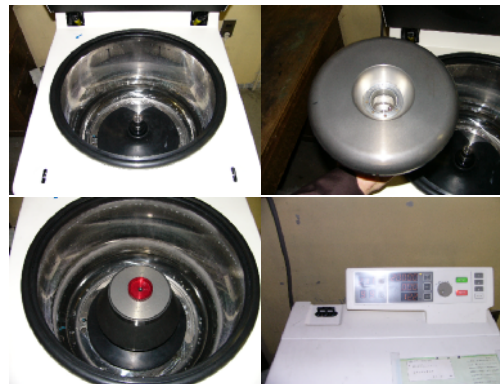

- (4) Set the tubes with the suspension (equally weighed) diagonally into the rotary part.

- (5) **Check and confirm the rotary part and machine's cover is connect with O-ring**, Close the rotary part by its cover then close the machine's cover by pressing it from the both side's hooks and set the working conditions, e.g. rotation speed, time and temperature. **Press start.**

- (6) At the finishing, take out your samples, clean and wipe out the inside wall and the rotary part then cover the rotary part and close the machine. Turn off the switch of the machine and of the main board. Remove the sample from the tubes and clean it nicely and keep those in the proper place. **Please don't forget to write your notes and remarks.** Keep the balance in unbalanced way as shown below: